

Pipetting the **SECRET CODE**

	1	2	3	4	5	6	7	8	9	10	11	12
A												
B												
C												
D												
E												
F												
G												
H												

Pipette the following amounts in the designated area within the 96 well plate. All values given below are in microliters (µl). Record a diagram of your plate using colored pencils in the chart above. Be sure to show your plate to an instructor once you are finished to record your finish order and scores. *Secret Code* _____ *Score* _____ *Accuracy Score* _____

Cell	Red	Blue	Green
D4	0	75	175
E4	0	85	165
F4	0	95	155
C8	140	110	0
F8	150	100	0
A12	55	0	195
G12	57	0	193
B11	62	0	188
C11	71	0	179
D11	91	0	159
E11	99	0	141
F11	101	0	149
C10	111	0	139
E10	121	0	129

Cell	Red	Blue	Green
C7	50	200	0
D7	60	190	0
F7	70	180	0
C6	80	170	0
E6	90	160	0
F6	100	150	0
C5	110	140	0
F5	120	130	0
C3	0	50	200
F3	0	60	190
C2	0	70	180
F2	0	75	175
D1	0	80	170
E1	0	85	165

D9	125	0	125
-----------	-----	---	-----

F1	0	90	160
-----------	---	----	-----